



FACSDiva Levey-Jennings Importer Instructions

1. Open the LJ Importer (.xlsm)


Double-click the **LJ Importer** file. Click the **Enable Content** button if the Security Warning below presents itself.

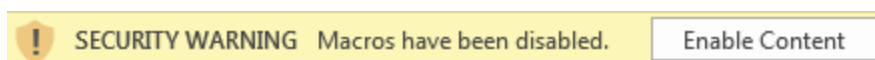

When opening the **LJ Importer** file from a server, the Security Warning below will be displayed. Note that answering Yes or No does not affect the following steps.

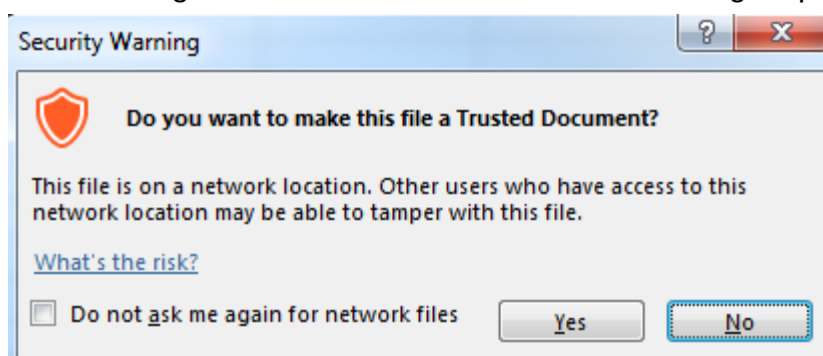


2. LJ Importer Buttons

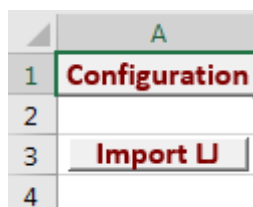

a) Configuration Button

Click on the **Configuration** button to select the Configuration.cst file those data you would like to load, then click the Open button. Alternatively, you can double-click on the Configuration.cst file to bypass clicking on the Open button.

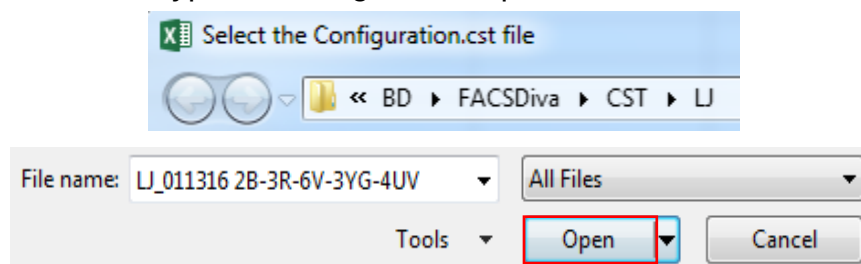


WORK-FLOW

Specialized Flow Cytometry Solutions

The configuration name will appear in cell **A2**.

	A	
1	Configuration	
2	2B-3R-6V-3YG-4UV.cst	
3	Import LJ	

Note that if the Configuration.cst file name is manually modified or if the file is moved or deleted after being selected, the following error message will be displayed:

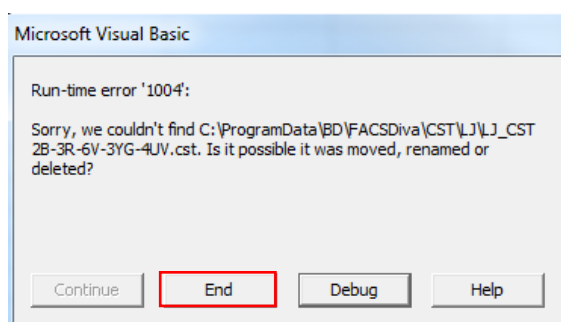


Click End to clear the error.

b) Import LJ Button

Once the Configuration.cst file has been selected, click the **Import LJ** button to transfer and parse the data to the **Data tab** (1st tab).

	A	B	C	D	E	F	
1	Configuration	Date_Time	Cytometer_Name	Cytometer_Type	Cytometer_SerialNo	Cytometer_Input_Device	Bead
2	eneron 2B-3R-6V	2015-11-24T09:03:14.9279257-08:00	LSRFortessa	LSRFortessa	1	Manual	CST Se
3	Import LJ	2015-12-18T14:58:27.9906291-05:00	LSRFortessa	LSRFortessa	1	HTS	CST Se
4		2016-01-06T15:15:30.4300010-05:00	LSRFortessa	LSRFortessa	1	HTS	CST Se

The graphs located in the **Charts tab** (2nd tab) will automatically populate. These graphs serve as examples, change them to fit your preferences.

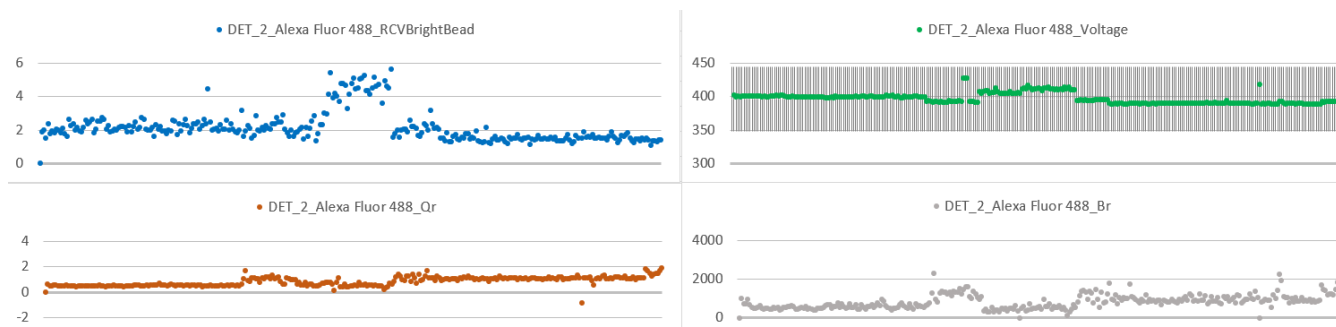

WORK-FLOW


Specialized Flow Cytometry Solutions

3. Modifying the Macro

The **LJ Importer** file contains 2 macros:

- Import_LJ_Click()
- ConfigFetch_Click()

To view and modify the code of these macros, go to the Developer tab and click View Code.

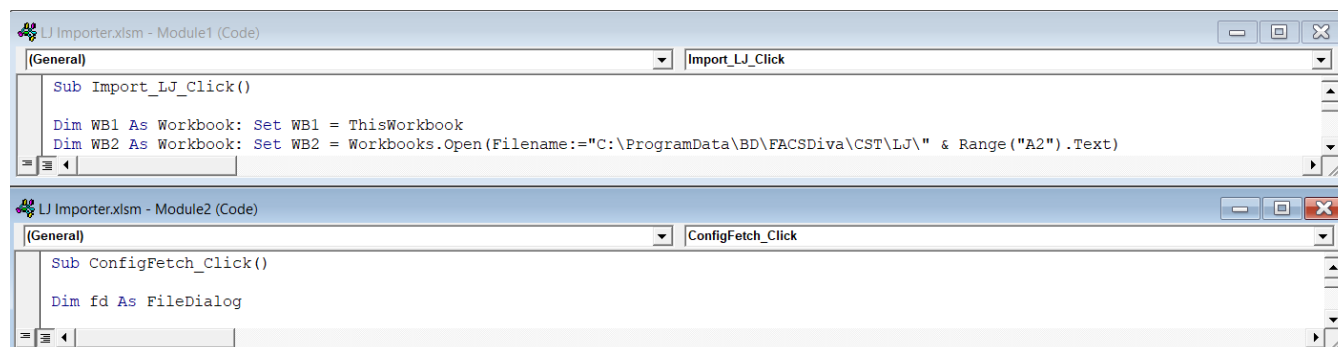

a) Checking data offline

The LJ CST files are located at: **C:\ProgramData\BD\FACSDiva\CST\LJ**

Several options are available to work with the data away from an instrument:

- Transfer the LJ CST file to a server
- Access the LJ CST file remotely
- Load the LJ CST file in Excel at the instrument, save the document then move it to an offline computer...

b) Testing the LJ Importer on a computer with no FACSDiva installation.

Create the following path: C:\ProgramData\BD\FACSDiva\CST\LJ

Place the **LJ Test** file provided in the zip package in the LJ folder.