

WORK-FLOW


Specialized Flow Cytometry Solutions

ForeCyt Data Store Script Instructions

1. The IntelliCyt Service

The **IntelliCyt Service** is set to automatically start when Windows boots up and is necessary to the proper functioning of the ForeCyt software.


However, occasionally, the service does not start or stops unexpectedly.

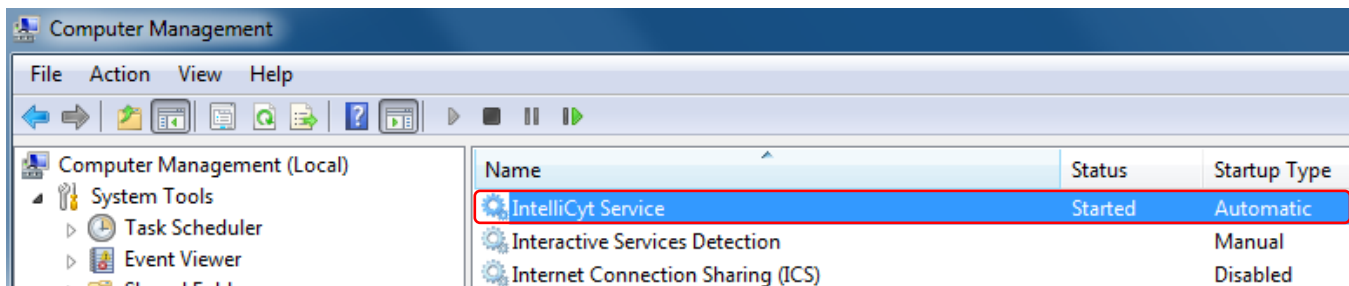

2. The Error

If the **IntelliCyt Service** is stopped, the following error will occur when launching the ForeCyt software.

Name	Status	Startup Type
IntelliCyt Service		Automatic

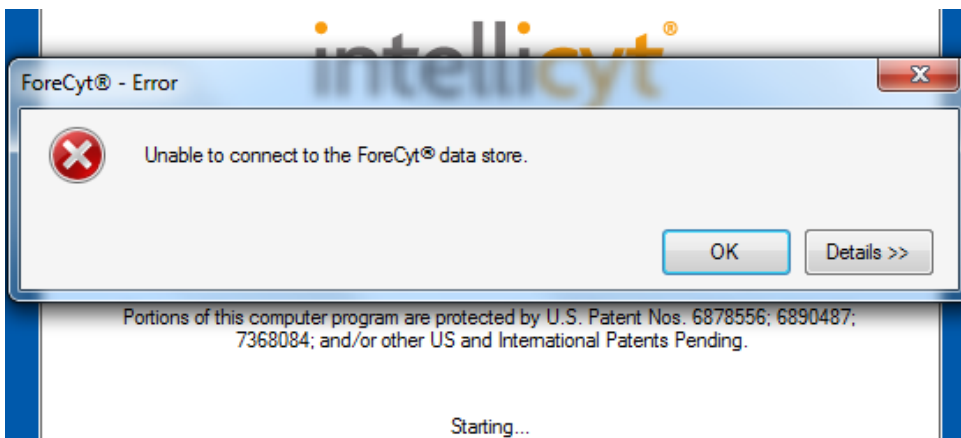

Error Detail:

Could not connect to net.tcp://localhost:8001/Integrated. The connection attempt lasted for a time span of 00:00:02. TCP error code 10061: No connection could be made because the target machine actively refused it 127.0.0.1:8001.

3. The Fix (requires administrator privileges)

Right-click the **ForeCyt Data Store Fix script** (.bat file) and select Run as administrator. The script simply starts the **IntelliCyt Service**.

WORK-*FLOW*

Specialized Flow Cytometry Solutions

Note: on an offline analysis PC if the **IntelliCyt Service** is stopped, the error appears as follows:

